# MicroTally™ Swab


*Improve Sampling Improve Quality*

Step 2

Step 3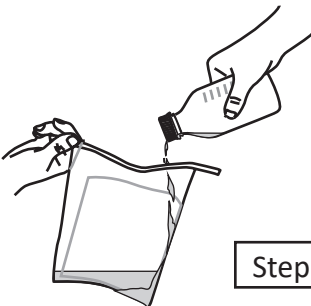

Step 6

1. Personnel responsible for collecting samples must wear plastic sleeves and gloves (not included). Prior to collecting each sample, personnel must sanitize sleeves and gloves with an approved sanitizer. Once gloves are sanitized, care should be taken so that gloved hands do not contact anything but the MICROTALLY™ Swab and product being tested.
2. Sampling can begin once the desired quantity of material is achieved within the combo. Aseptically remove the Manual Sampling Device (MSD) MICROTALLY™ Swab from the sample bag, and swab the entire top surface of the combo. Make sure to use enough pressure to ensure that any bacteria present are dislodged from the product and captured within the MICROTALLY™ Swab.
3. Using one side of the MSD MICROTALLY™ Swab, swab one half of the meat exposed on the top surface of the combo, flip the MICROTALLY™ Swab and swab the remaining half of the top surface of the combo. Sample collection is conducted for approximately 1.5 minutes.
4. After sample collection, place the used MICROTALLY™ Swab, with an additional horizontal fold, into original bag. Label bag following company procedure.
5. Repeat steps as necessary to sample additional combos; each combo will require a new MICROTALLY™ Swab. To prevent cross-contamination, gloves should be changed after each combo.
6. To analyze the samples, enrich the MICROTALLY™ Swabs with 200 mL of appropriate nutrient broth and analyze via the same procedures normally utilized for combo lot testing.

## FREMONTA CORPORATION

FOR TECHNICAL SUPPORT OR TO PLACE AN ORDER

VISIT [WWW.FREMONTA.COM](http://WWW.FREMONTA.COM) OR EMAIL [ORDER@FREMONTA.COM](mailto:ORDER@FREMONTA.COM)

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